


Cooling Incubator

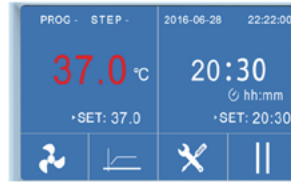


Advanced Color Intelligent Touch Screen Control

- 5.0 Inch Touch screen, intelligent touch control, real-time display of parameters, simple and convenient operation.
- Quick setting of temperature, time and other parameters can be carried out. It has the function of program setting.
- It can be programmed in 7 segments, 63 steps, 9 steps and 1-99 hours and 59 minutes per segment.
- With touch screen automatic locking function




Professional LCD Controller (L Series)

- Intelligent program control with high brightness LCD screen, button operation and real-time display of parameters.
- Quick setting of temperature, time and other parameters can be carried out.
- It has the function of program setting. It can be programmed in 7 segments, 63 steps, 9 steps and 1-99 hours and 59 minutes per segment.

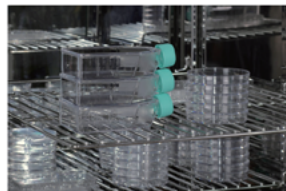

Anti-skid and anti-overturning shelf design

- Shelf has anti-inclination function to avoid falling off when shelf is pulled out.
- Shelf has anti-lock function, the shelf will be locked when it is pulled out half, so as to avoid the loss caused by the direct slide of the shelf.

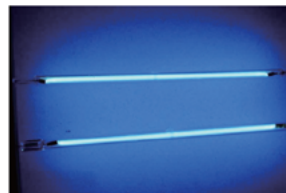

Ultraviolet sterilization system (option)

- Ultraviolet lamp is located on the upper wall of the box body. It can disinfect the box body regularly and kill the plankton bacteria in the circulating air of the box body effectively so as to prevent the pollution during cell culture.



Professional duct design

- With a unique duct design, the use of large impeller fans, fast temperature rise and fall, stable temperature control, good temperature uniformity.
- The circulating fan has adjustable speed, middle and low grades, so as to avoid the volatilization of samples due to excessive air flow during the test.
- Famous brand fan, high efficiency, low noise, energy saving and environmental protection.



Stainless steel inner liner

- 304 stainless steel, mirror polishing, strong anti-corruption ability.
- Large arc angle design, no dead angle, easy cleaning and maintenance.

Refrigeration system

- International brand compressors, with high refrigeration efficiency and low noise, shorten cooling time by 40% compared with traditional cryogenic equipment, save energy and protect environment.
- Minimum temperature up to -10°C, with refrigeration and heating functions.
- Automatic switching of refrigeration and heating, high control accuracy and fast temperature stability.

Reserved test hole

- External detection sensor can real-time detect the temperature inside the box through the test hole to ensure that the temperature inside the box is accurate and reliable.

Cooling Incubator

Humanized design

- With casters, flexible and convenient to move.
- With reservation and timing functions, there is no need to wait, which effectively improves the efficiency of the experiment.
- Available in Chinese and English menus to meet different language needs, with °C / °F conversion.

Safety


- Protection on instruments: Comply international standard secondary temp limiting alarm system. When the heating is out of control or exceeds the maximum limit temperature, the heating is automatically cut off with sound and light alarm, ensure operating is safe without any accident.
- Protection on key components: Key components are equipped with over-current, over-temperature, overload and other safety protection to prevent accident.
- Protection on samples: When the temperature inside the box is higher or lower than the set temperature, the alarm will start to cut off the heater, with sound and light alarm to remind the operator to protect the sample from normal test without accident.
- Protection on operator: The cabinet and door are specially insulated to make the surface temperature of the cabinet low, ensuring the operator's safety and no accidents.
- Breakdown message provided: When the instrument breakdown, the message will shown on the screen to help operator clearly check.

Convenient data processing

The touch screen type standard with a USB interface, which can record the change status of the temperature parameters. The LCD screen is option for USB interface.

- It can be equipped with RS232 data interface, which can realize remote control of the machine through software (option).
- RS-232 and USB can be selected one of them.

Cooling incubator Structure diagram

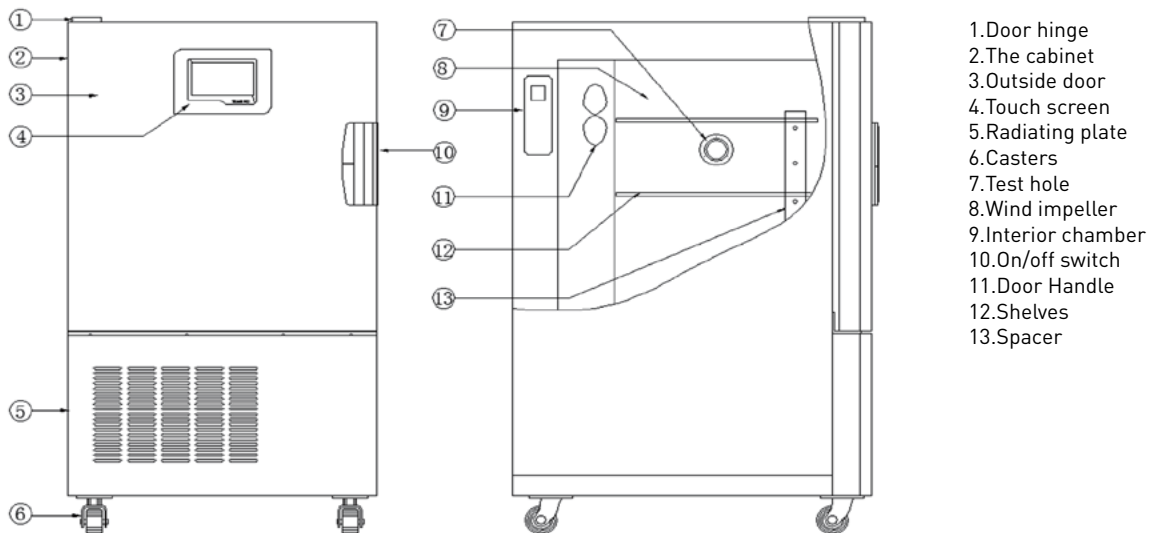

Cooling incubator



Model	BC-60 BC-60L	BJ-60 BJ-60L	BC-120 BC-120L	BJ-120 BJ-120L	BC-250 BC-250L	BJ-250 BJ-250L	BC-500 BC-500L	BJ-500 BJ-500L
Capacity (L)	68L		120L		252L		495L	
Temp range	-10~80°C							
Display Resolution	0.1°C							
Uniformity	±1.0°C							
Timing Range	1~5999mins							
Max number of shelves	2(10)		3(14)		3(16)		4(16)	
Max load per shelf	20Kg							
Net Weight	85Kg		100Kg		120Kg		220Kg	
Interior Dimension W×D×H(mm)	400×380×450		500×400×600		550×600×750		670×725×1020	
Exterior Dimension W×D×H(mm)	545×670×1000		645×690×1150		695×890×1300		850×1095×1935	
Exterior Dimension W×D×H(mm)	545×670×1000		645×690×1150		695×890×1300		850×1095×1935	
Power Consumption	AC220V/50Hz							
UV lamp	Optional	Standard	Optional	Standard	Optional	Standard	Optional	Standard

BC/BJ series touch screen controller
BC series Cooling incubator

BC/BJ-L series LCD controller
BJ series Cooling incubator with uv lamp

Heating incubator




Advanced Color Intelligent Touch Screen Control

- 4.3 Inch Touch screen, intelligent touch control, real-time display of parameters, simple and convenient operation.
- Quick setting of temperature, time and other parameters can be carried out.
- It has the function of program setting. It can be programmed in 7 segments, 63 steps, 9 steps and 1-99 hours and 59 minutes per segment.
- With touch screen automatic locking function to prevent non-related personnel from disoperation.

Professional LCD Controller (L Series)

- Intelligent program control with high brightness LCD screen, button operation and real-time display of parameters.
- Quick setting of temperature, time and other parameters can be carried out.
- It has the function of program setting. It can be programmed in 7 segments, 63 steps, 9 steps and 1-99 hours and 59 minutes per segment.

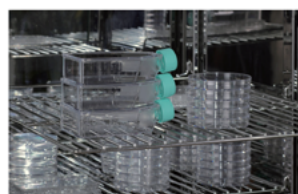


Professional anti-skid and anti-overturning shelf design

- Shelf has anti-inclination function to avoid falling off when shelf is pulled out.
- Shelf has anti-lock function, the shelf will be locked when it is pulled out half, so as to avoid the loss caused by the direct slide of the shelf.



Temperature test hole

- External detection sensor can real-time detect the temperature inside the box through the test hole to ensure that the temperature inside the box is accurate and reliable.

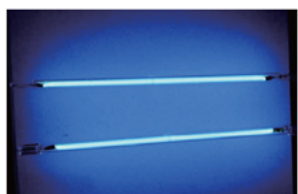

Ultraviolet sterilization system (option)

- Ultraviolet lamp is located on the upper wall of the box body. It can disinfect the box body regularly and kill the plankton bacteria in the circulating air of the box body effectively so as to prevent the pollution during cell culture.



Temperature limit alarm system

- When the temperature control system is out of control and the temperature exceeds the limit temperature, the heating is automatically interrupted to ensure the safety of testing.



Stainless steel inner liner

- 304 stainless steel, mirror polishing, strong anti-corruption ability.
- Large arc angle design, no dead angle, easy cleaning and maintenance.



Professional duct design

- With a unique duct design, the use of large impeller fans, fast temperature rise and fall, stable temperature control, good temperature uniformity.
- The circulating fan has adjustable speed, middle and low grades, so as to avoid the volatilization of samples due to excessive air flow during the test.
- Famous brand fan, high efficiency, low noise, energy saving and environmental protection.



Sample temperature sensor (optional)

- Optional sample temperature sensor, real-time detection of sample temperature, so that the temperature is more real and reliable.
- Double Convection Model
- Fan power can be adjusted from 0% (equivalent to natural convection) - 100% four gears, which can meet the training requirements of different airflow.

Heating incubator

Humanized design

- Adopting direct heating system, the heating speed is fast and the control precision is high.
- PID control technology, high control precision and small temperature fluctuation.
- Available in Chinese and English menus to meet different language needs, with °C / °F conversion.
- The height of the shelf is adjustable to meet different cultivation needs.
- Professional stacking foot design, the machine can be stacked, saving lab space and improving the use efficiency (same specification).

Convenient data processing

- The touch screen type standard with a USB interface, which can record the change status of the temperature parameters. The LCD screen is option for USB interface.
- It can be equipped with RS232 data interface, which can realize remote control of the machine through software (option).
- RS-232 and USB can be selected one of them.

Safety

- Protection on instruments: Comply international standard secondary temp limiting alarm system. When the heating is out of control or exceeds the maximum limit temperature, the heating is automatically cut off with sound and light alarm, ensure operating is safe without any accident.
- Protection on key components: Key components are equipped with over-current, over-temperature, overload and other safety protection to prevent accident.
- Protection on samples: When the temperature inside the box is higher or lower than the set temperature, the alarm will start to cut off the heater, with sound and light alarm to remind the operator to protect the sample from normal test without accident.
- Protection on operator: The cabinet and door are specially insulated to make the surface temperature of the cabinet low, ensuring the operator's safety and no accidents.
- Breakdown message provided: When the instrument breakdown, the message will shown on the screen to help operator clearly check.

Model	BI-35T BI-35TL	BI-55T BI-55TL	BI-120T BI-120TL	BI-200T BI-200TL
Capacity (L)	34L	54L	124L	222L
Temp range	Amb +5~80°C			
Display Resolution	±0.2°C			
Uniformity(at 37°C)	±0.8°C			
Display Resolution	0.1°C			
Max number of shelves	2(5)	2(9)	2(12)	2(16)
Max load per shelf	20KG			
Net Weight	44Kg	53Kg	79Kg	108Kg
Timing Range	1~5999mins			
Interior Dimension W×D×H(mm)	320×300×320	400×330×380	520×450×490	650×500×650
Exterior Dimension W×D×H(mm)	610×520×580	690×468×640	810×588×755	940×658×910
Electrical Requirement	AC220V,50Hz			
Power Consumption	250W	300W	550W	700W

Heating incubator

Technical Parameters (Air Forced)

Model	BI-35F BI-35FL	BI-55F BI-55FL	BI-120F BI-120FL	BI-200F BI-200FL	BI-400F BI-400FL
Capacity (L)	35L	67L	146L	253L	420L
Temp range	Amb +5~80°C				
Display Resolution	±0.2°C				
Uniformity(at 37°C)	±0.8°C				
Display Resolution	0.1°C				
Max number of shelves	2(5)	2(6)	2(10)	2(16)	23
Max load per shelf	20Kg				
Net Weight	48Kg	56Kg	82Kg	119Kg	160Kg
Timing Range	1~5999mins				
Interior Dimension W×D×H(mm)	320×300×320	400×305×415	520×430×530	650×500×650	1000×800×530
Exterior Dimension W×D×H(mm)	610×550×545	690×560×640	810×685×755	940×755×875	1285×980×750
Electrical Requirement	AC220V,50Hz				
Power Consumption	300W	350W	600W	700W	1000W

BI-F series touch screen controller

BI-FL series LCD controller

- The outer dimensions do not include hinges, handles, gauges, wires, etc., and the internal dimensions do not include the duct section.
- All test data are tested in a 25°C environment.
- If the maximum temperature is to reach 300°C , please note in advance before ordering.